

# bcAnalyze - Manual

Version: August 2016, Version 3.0

Author: Dr. Volker Runkel

## Contents

<b>1. Introduction</b>	<b>3</b>
<b>2. Installation</b>	<b>3</b>
<b>3. Quick start</b>	<b>3</b>
<b>4. Basics</b>	<b>5</b>
<i>4.1. Sound definitions</i>	<i>5</i>
<i>Sound</i>	<i>5</i>
<i>Frequency</i>	<i>5</i>
<i>Microphone</i>	<i>5</i>
<i>Sample rate</i>	<i>5</i>
<i>Amplitude resolution</i>	<i>6</i>
<i>Sound Pressure Level, override (Clipping)</i>	<i>6</i>
<i>Frequency analysis</i>	<i>6</i>
<b>5. First steps</b>	<b>7</b>
<i>5.1. Open files</i>	<i>7</i>
<i>5.2. Zooming and selecting sound</i>	<i>7</i>
<b>6. Analyzing signals</b>	<b>8</b>
<i>6.1. Spectrum</i>	<i>8</i>
<i>6.2. Sonogram</i>	<i>9</i>
<i>6.3. Call overlay</i>	<i>11</i>
<i>6.4. Waveform statistics</i>	<i>12</i>
<b>7. Finding and measuring calls automatically</b>	<b>12</b>
<i>7.1. The call finder and its requirements</i>	<i>12</i>
<i>7.2. Working with the automatic call finder</i>	<i>13</i>
<b>8. Play calls</b>	<b>13</b>
<i>8.1. Time-delay player</i>	<i>14</i>
<i>8.2. Virtual heterodyne detector</i>	<i>14</i>
<b>9. Sound manipulation</b>	<b>14</b>
<i>9.1. Changing samplerate</i>	<i>14</i>
<i>9.2. Amplify</i>	<i>14</i>
<i>9.3. Generate</i>	<i>14</i>
<i>9.3.1. Silence, white noise, sinus</i>	<i>14</i>
<i>9.3.2. Generate bat call</i>	<i>14</i>
<i>9.4. Filter</i>	<i>15</i>
<b>10. bcCalls documents</b>	<b>15</b>
<b>11. CSV documents</b>	<b>16</b>

<b>12. Preferences</b>	<b>16</b>
<b>12.1. General</b>	<b>16</b>
<b>12.2. Display</b>	<b>16</b>
<b>12.3. Call finder</b>	<b>16</b>
<b>12.4. Sonogram</b>	<b>16</b>

# 1. Introduction


With bcAnalyze you have one of the most modern tools for analyzing bat echolocation and social calls. While the application was fine-tuned for using bat corder recordings other files in wave-format can be read and analyzed nearly as well. The application focuses on a fast display and analysis of sound files embedded in an easy to use user interface. A special algorithm for finding bat calls within recordings allows the automatic extraction of call parameters. The following pages cover bcAnalyze 3 Light as well as the Pro version. Pro chapters and functions are marked as such.

## 2. Installation

bcAnalyze requires an Apple Mac running OS X 10.9 or newer. We recommend memory of at least 4GB. Both versions, light and pro, are available through the Mac App Store and installed from there automatically after purchase.

## 3. Quick start

If you have worked with other sound analysis tools before, you recognize bcAnalyze is built in a similar way. After opening a sound file (batcorder raw or windows wave), you see an waveform or oscillogram display of the sound. Above the wave form a sonogram for quick call display is shown. If the sound file was already analyzed with the call finder (for example in bcAdmin) and a bcCalls file was stored, the found calls are displayed as well as overlay over the sonogram (can be disabled).

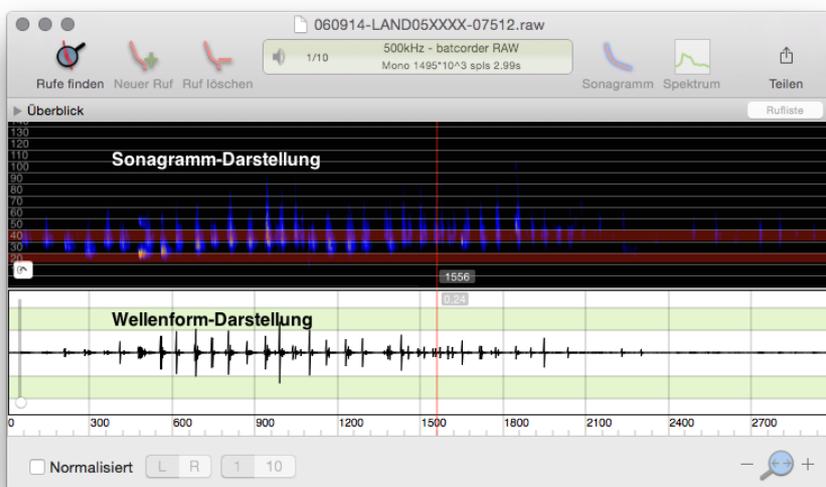


After selecting a piece of sound you can calculate and show a spectrum display or a measurement sonogram. If found calls are available the tab key quickly browses through the call list while ⌘+R creates a sonogram of the selected call. See more on the features of bcAnalyze on the following pages.

## 4. Basics

### 4.1. Sound definitions

#### Sound

Sound is an oscillation of air pressure which spreads as a wave with a velocity of ca. 340 m/s. An oscillation of one specific frequency is called a pure tone, which results in a sinus wave when outlined over time. Such an illustration is called wave display or oscillogram.

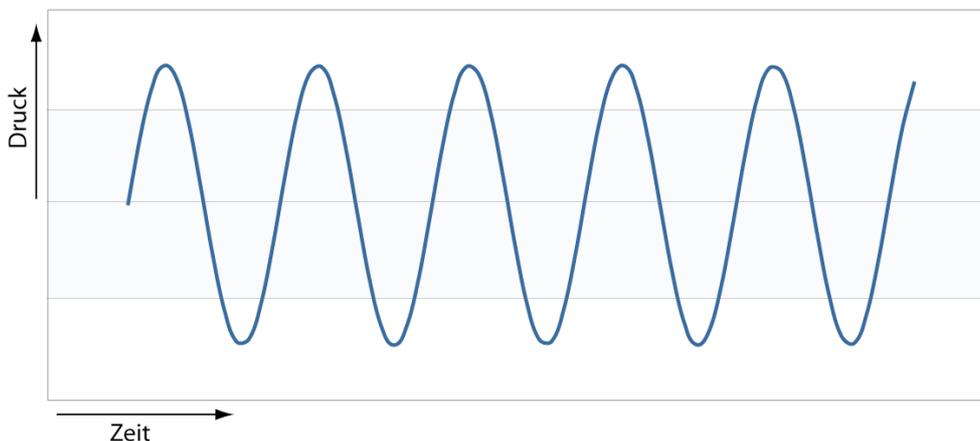

#### Frequency

The frequency of sound is defined by the number of oscillations per second [Hz - Hertz]. Depending on the frequency sound can be classified as infrasonic sound (below 16 Hz), audible sound (16 Hz up to 20 kHz, audible for humans), and ultrasonic sound (20 kHz up to 1 GHz). In contrast to a pure tone, sound (in terms of noise) is composed of many frequencies that change over time. A tone is a periodic signal generated by several pure tones.

#### Microphone

A pressure sensor (microphone) detects changes in air pressure and transforms these into electrical pulses. These pulses ideally (exact linear transformation) form a wave as explained above. This is still an analogue signal (continuous signal), which has to be transcribed in discrete time and amplitude steps in order to digitalize it.

#### Sample rate

The sample rate defines the time interval for reading out the signal. The sample rate has to be significantly higher than the frequency of the respective signal. For an accurate digital reproduction of an acoustic signal, the sample rate has to be twice as high as the maximal frequency of the analogous signal (Nyquist-Shannon Sample Rate Theorem). An audio CD for example is recorded with a sample rate of 44.1 kHz in order to reproduce human audible frequencies from 50 Hz up to 20 kHz. For bat calls the sample rate should be at least

250 kHz because echolocation calls can reach up to 125 kHz. We recommend a sample rate of 500 kHz for the automated analysis with bcAnalyze.

### **Amplitude resolution**

This attribute denotes the possible amount of amplitude values for digitalization. For a resolution of  $N$  exactly  $2N$  values have to be available, e.g. resulting in 65536 steps for 16 Bit. Generally, in acoustic digitalization each sample becomes digitalized with at least 16 Bit resolution, which corresponds to a theoretical dynamic of 96 dB. We recommend to apply at least 16 Bit for a feasible analysis of ultrasonic signals.


### **Sound Pressure Level, override (Clipping)**

The sound pressure level (loudness) can be specified linearly (pressure in Pascal or voltage in Volt). However, it is more common in acoustics to use a logarithmic scale: Decibel dB, which describes the loudness with the common logarithm of the ratio of two levels. Therefore it is a relative unit without dimension, representing an increase with positive and an attenuation with negative values. dB specifications are often related to a standardized reference value: dB SPL is referenced to the human auditory threshold at 1 kHz and indicates how much louder a signal is as compared to this value. The great range between quiet and loud can be comfortably expressed by this unit. For instance, a difference of 80 dB between two signals relates to a voltage ratio of 1:10.000. As a convenient rule of thumb: a doubling of the sound pressure level (respectively voltage) results in 6 dB, a ten-fold increase in 20 dB difference.

An override results if the analogue signal is louder than the input of the digitalization device (AD-Converter). The sinus wave is clipped and the signal gets adulterated. This also applies to analogue amplifiers as they cannot produce a linear amplification (distortion). In frequency analysis this causes deceptive frequencies.

### **Frequency analysis**

There are several methods for an analysis of sound signals. A simple and feasible method is the zero crossing analysis: it calculates the oscillations per second (frequency) by counting how often the zero point is crossed and establishes a periodogram. With many consecutive measures a frequency progression over time can be constructed (see Figure to your right-hand side). The Fourier Transformation is the most common method to analyze

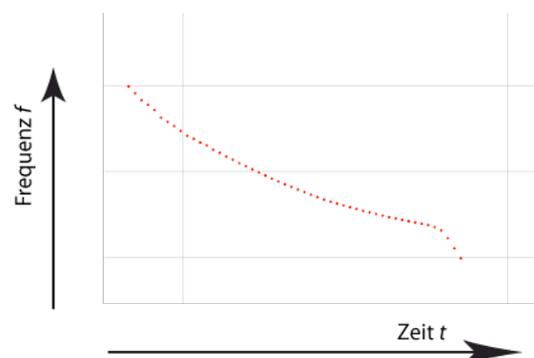

an acoustic signal which is composed of several frequencies. This elaborate technique separates a periodic signal to its single frequency components. bcAnalyze applies a slightly

less complex method, the Fast Fourier Transformation (FFT) for the calculation of spectra and sonograms. Compared to a spectrum, which illustrates the frequency classes of a signal, a sonogram also depicts the temporal pattern of these frequencies.

## 5. First steps

### 5.1. Open files

Using the Open command or the Recent Documents menu, both located in the File menu you can open files from hard disk. Alternatively bcAnalyze supports drag and drop from the Finder. If bcAnalyze is set as standard application for a file type, it opens upon double-clicking a sound file in a Finder window.

Currently support for bat corder .raw and wave .wav files exists. Note that compressed wave formats are not supported. If a wave file contains two channels (stereo), only one is displayed at a time. You can easily switch between left and right channel using the channel chooser below the oscillogram display. Wave files with a sample rate below 196 kHz are interpreted as time-delayed recordings automatically.

After opening the file the waveform shows a linear display of the recorded sound wave. Signals that reach the upper or lower boundary of the display are distorted. Below the waveform display to the left a checkbox allows to Normalize sound display based on the visible samples. In addition to normalizing the sound one can also use the slider control to change the y-axis magnification. The x-axis shows the time and is labeled in milliseconds. If call were found or exist for a recording bars on the x-axis indicate their position. Above the waveform a sonogram display is visible.

**Pro:** To save changes to a file you can use File - Save as... and choose to store in raw or wave format.

### 5.2. Zooming and selecting sound

Per default the full sound is displayed as waveform and sonogram directly after loading the sound file. To zoom into the sound to get a more detailed view you can either use the +/- keys or the buttons adjacent to the small magnifying glass symbol to the windows lower right. A click on the magnifying glass zooms out and shows the full sound again. If you hold the shift key while clicking the magnifying glass the current selection is zoomed to.

When moving the cursor over the waveform the current time and voltage value is displayed. With the left mouse key held down a selection can be drawn in the waveform, indicated by the red rectangle. A selection will be used to calculate spectrum, sonogram or do a manipulation of that range of sound samples. The duration of the current selection is shown rounded to milliseconds. Using the command Set selection from within the Edit menu allows a sample exact selection of sound. The Edit menu features a deselect and a select all

command as well. An alternative is to create the selection by drawing a rectangle within the sonogram. This displays not only the duration of the selection but also lowest and highest frequency of the selected range.


The sonogram display as such can be adjusted quickly by moving the mouse over the arrow-up button and choosing either a slider for color changes or one of the two options. The one called Calls toggles display of call measurements. These are, if activated, plotted above the sonogram as an overlay. The second toggles display of the sonogram as such. The sonogram features two bands helping quickly identifying frequency ranges of recorded calls. These bands can be adjusted in the application preferences. By pressing the shift key a help line is added at the current mouse position. Another press of shift disables the line again.

## 6. Analyzing signals


### 6.1. Spectrum

A power spectrum is calculated using a fast fourier transformation and shows the amplitude or power distribution of the frequencies contained within the analyzed signal. A spectrum doesn't contain any temporal information. Usually it is calculated over a single analysis window defined by the sound selection. The spectrum, if used correctly, is rather accurate regarding frequency measurements.

The window function defines a weighting of samples within the analysed block. While a rectangle window weights all samples with the same, full value, the other window functions have a bell shaped distribution of weights. Thus, parts of the samples located at the beginning and end of the window are less pronounced than the middle part. The larger the window size the higher is the frequency resolution. bcAnalyze displays the resolution when calculating spectra. Spectrum calculation is working for sample windows of a size of  $2^N$  samples. bcAnalyze uses zero-padding to fit the selected sound into such a window. So first you have to select the sound range that you want to calculate a spectrum from. Usually that will be a single bat call. Choose Spectrum either from the toolbar or from the Analyse menu and the spectrum will be displayed in a new window. The graph shows a logarithmic display of power and a li-

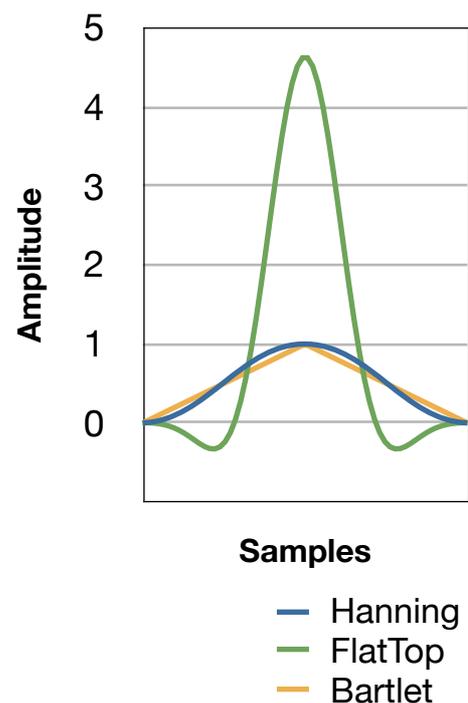


near display of frequencies. You can choose the currently used window function from the

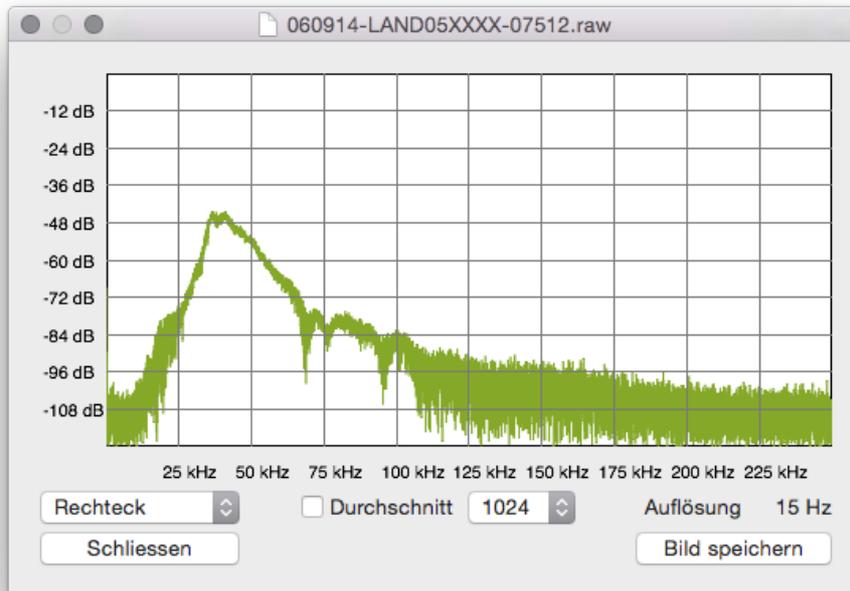


popup menu. Moving the mouse over the spectrum activates a cursor that displays frequency and amplitude/power. For a smoother display you can use a smoothing function that averages based on your selection.

## 6.2. Sonogram


A sonogram is generated by calculating many spectra over very short sample ranges and combine them into a quasi two-dimensional image showing frequency over time. The frequencies amplitude is color-coded adding a third dimension. In contrast to a spectrum display the optimal window function is not a rectangle but one that has either a good temporal or frequency response, or both. These types reduce the sidelobes (picket fences). The available window types in bcAnalyze have all pros and cons regarding the resolution, the best window function is the 7th term harris window. The following images show three different window types for the same signal.

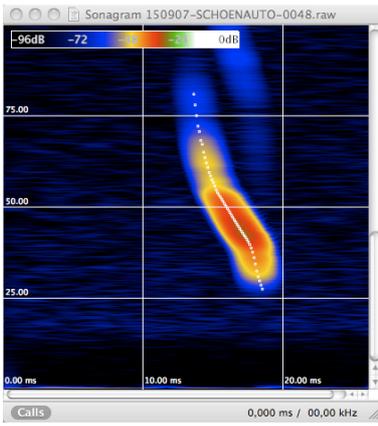

Hanning

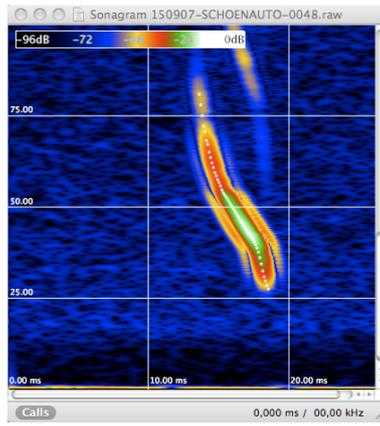


FlatTop

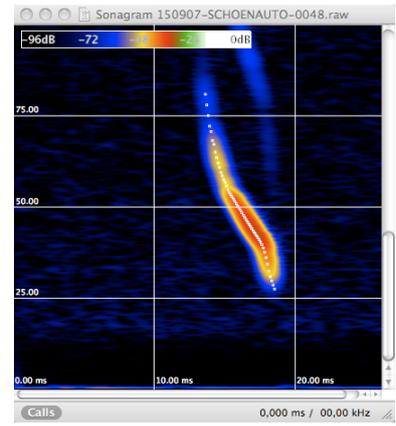

7th term Harris

Frequency and temporal resolution of a sonogram is influenced by various parameters. Mainly the FFT window size and the overlap both have the highest impact. The FFT size is the number of samples analyzed per single calculation. The higher its value, the finer the frequency resolution, while temporal resolution decreases. The overlap indicates how many samples each window start are spaced. High values produce high temporal resolutions. Whenever a sonogram is opened in an extra window in bcAnalyze the resolution is displayed shortly. Apart from window size and type the sonogram display can be adapted by using the color popup in the upper right as well as the sliders.

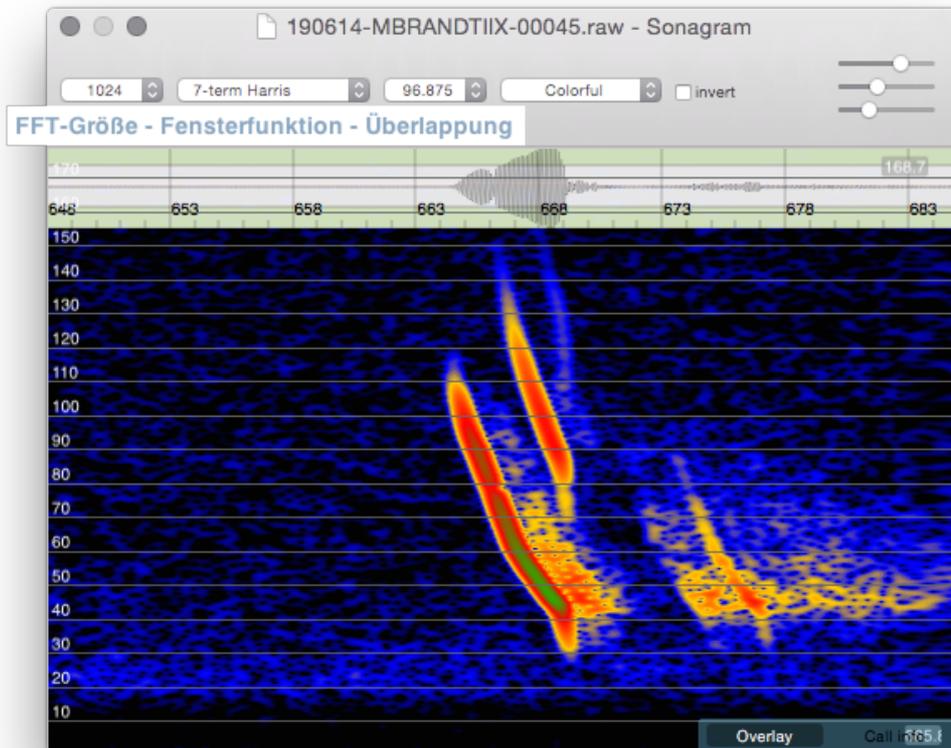


Moving the mouse over the sonogram show the current frequency and time at the mouse cursor. While holding down the left mouse key one can draw a selection. The selection features duration as well as upper and lower frequency. This helps to quickly get start-, end-frequency and duration of a bat call.

The button control in the lower right toggle two kind of measurement overlays if measurements are available for the visible call.

If you have changed sonogram settings you can store them as default by doing a right click on the sonogram and choosing the appropriate item from the context menu. The same menu offers various other possibilities like saving the sonogram to an image or an html page. The latter requires you to choose a folder to which an index file and the image is saved. If the folder already contains an export produced by bcAnalyze the new sonogram is attached to the existing page. That way you can easily build a single web page featuring a multitude of sonograms.

**Pro:** The pro version offers the possibility to copy a selected sonogram area and paste it into another sonogram for comparing calls. The image is also available in the general paste-board for other applications. After pasting such a sonogram piece you can move it by either using the mouse, holding the shift key locks movement to one axis, or by the arrow keys.

The arrow keys can be used together with shift to increase the movement steps. Using page up or down increases or decreases the opacity of the pasted sound piece. The Delete backwards key deletes the pasted image again.

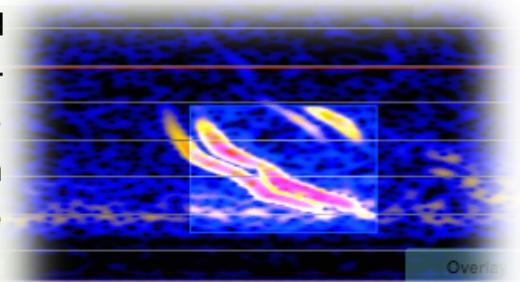

### 6.3. Call overlay

Another tool to help identifying bat species is an overlay of call measurements. For this all found calls are drawn into a rough grid in a way that the call ends are aligned. In addition you can overlay stored references as well as self generated references from your own call recordings using the bcCalls files.

**Pro:** You can open bcCalls files as own documents and copy & paste call data into one large reference bcCalls file for the call overlay. The pro version offers to open bcCalls files as separate documents.



#### 6.4. Waveform statistics

**Pro:** Located in the Analysis menu is a simple tool for evaluating waveform statistics of the currently selected sound. It lists DC offset as well as minimum and maximum amplitude of the selection. The results are shown as percentage and as dB value relative to the maximum possible amplitude. The average RMS (root mean square) value is given in addition.

## 7. Finding and measuring calls automatically

bcAnalyze offers a unique algorithm for finding bat calls within recordings. Found calls are automatically measured for species identification and values are accessible by the user as well. The calls are modeled by getting the instantaneous frequency for every 0.1 ms. This data is displayed as overlay over the calls in the sonogram displays.

### 7.1. The call finder and its requirements

The call finder and measurement extractor is based on a complex form of zero-crossing analysis combined with statistical tests. In a first step possible calls are evaluated using a choosable amplitude threshold. In the next steps all found calls are evaluated using statistics. That way badly recorded calls or non-call sounds are mostly eliminated. For the valid calls parameters like start- and end frequency as well as the call overlay data are extracted.

Some requirements should be met by the recordings for a robust call extraction:

- the sound file must contain ultrasonic signals > 14 kHz
- the call must be tonal with a single accented main harmonic

- the call must have a duration of at least 1.5 ms
- overlapping signals (call and echo for example) can not be extracted properly
- the call amplitude must at least reach the threshold given in - xx dB relative to maximum amplitude
- the recording background noise must be low respectively the signal-to-noise ration must be good
- loud background signals must not mask the bat calls

Some of the call finder parameters can be adjusted. Within the application preferences you find an area for the call finder. It allows adjustment of the call finder threshold. The lower the value (note -36 dB is lower than -30 dB), the more calls can potentially be found. The quality setting is used for the statistical tests of the call finder. The higher the value the lower the quality can be to get a found call. A value of 20 usually is optimal for all call situations. If you lower the value towards zero fewer calls will be found. If you work with bat corder recordings you should adapt the call finder threshold to the bat corder threshold value for optimal performance. The selected threshold can be displayed in the waveform view. Choose View from the main menu and Show threshold within there. It is displayed as a horizontal line.

## 7.2. Working with the automatic call finder

A recording can be searched for calls using the Call finder toolbar action or Find calls from the Analysis menu. If recordings were analyzed using bcAdmin the calls are already stored per recording in the .bcCalls file and are automatically read from there. For example this is a pair of files to expected after bcAdmin analyzed the recording:

```
250508-ZWEIM2XXXX-0004.raw
250508-ZWEIM2XXXX-0004.bcCalls
```

The found calls are listed as a table in the call list popup. There you can read out values like start- and end frequency as well as user defined modal frequency (FMod), Frequency of last millisecond (Flastms) or Frequency at lowest slope (FRmin).

**Pro:** After calls were found these are not automatically saved to disk. Using the export option in the Calls menu (main menu) will save the found calls to a bcCalls file.

## 8. Play calls

**Pro:** Some users prefer to listen to recorded calls. The integrated sound player either replays a time-delayed version of the recording or a heterodyne version. You can start the player either via the speaker symbol in the toolbars info field or using ctrl+space.

### 8.1. Time-delay player

**Pro:** This mode reduces the samplerate by a factor of 10 without altering the sound. Thus, the originally ultrasonic signals are played at a tenth of their recorded frequency and thus audible to us.

### 8.2. Virtual heterodyne detector

**Pro:** The heterodyne mode simulates a handheld heterodyne bat detector. You can choose the frequency and listen to your recording just like in the field.

## 9. Sound manipulation

**Pro:** Various tools for sound manipulations exist in bcAnalyze Pro. These allow to filter sound, change the samplerate or generate sounds.

### 9.1. Changing samplerate

**Pro:** The Edit menu holds an item for changing the samplerate. This can be done with or without changing the samples. In the first case the sound is resampled to the new sampling rate. If you lower the rate, frequency above half the new rate will be gone in the new sound. If you deactivate resampling the frequencies will shift according to the difference of old and new samplerate.

### 9.2. Amplify

**Pro:** The command Edit-Amplify increases or decreases the signal amplitude according to the given value.

### 9.3. Generate


**Pro:** You can generate silence, white noise, sinus signals and bat calls by choosing the appropriate option from the Edit - Generate menu.

#### 9.3.1. Silence, white noise, sinus

**Pro:** These functions allow generation of the selected signal-type for either the selection duration or a given duration. The signal is inserted at the waveforms insertion point.

#### 9.3.2. Generate bat call

**Pro:** You can create a bat call like signal that can be used with a bat lure for example. The tool allows you to „draw“ a call that can be then generated as sound including amplitude variations. You either have to create a new 500 kHz sound file or open one with 500 kHz samplerate. Choose from the Edit- generate menu the Bat call generator and a window opens up for drawing the bat call.

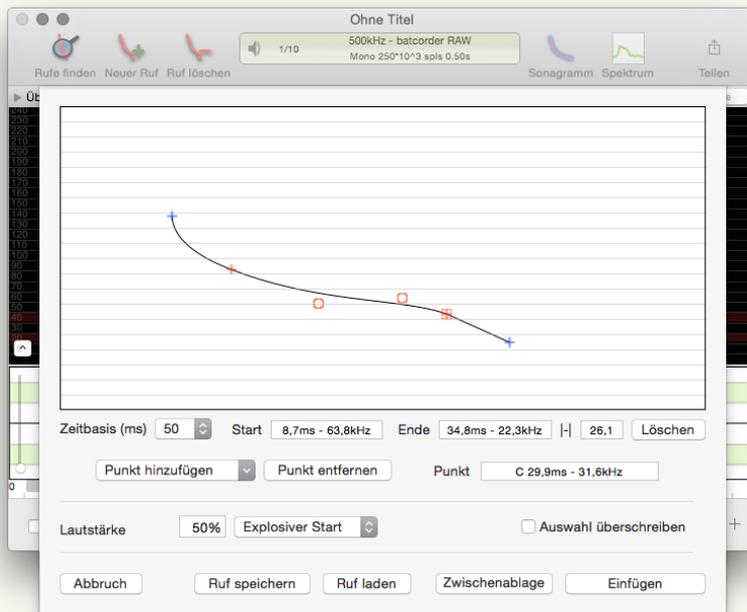

Designed calls can be saved to disk and reopened later again. Use the button Clipboard or Insert to generate the designed call.

#### 9.4. Filter

**Pro:** Simple biquad-filters allow bandpass-, low- and highpass filtering of selected sounds.

## 10. bcCalls documents

**Pro:** If calls are measured with bcAdmin the results are stored in .bcCalls files. Creating these from bcAnalyze is possible as well. These files can be opened just like a recording. The stored calls are listed in a table. Call overlays and optionally the measurements can be displayed for all calls. Calls can be copied between bcCalls documents. That way you can for example generate a file to be used with the call overlay feature of bcAnalyze.

#	Start	Dur	IPI	SF	EF	FMod	Flms	FRmin	Species	Prob
1	88,3	1,6	---	122,0	19,3	105	64,0	106,4	Mnat	100
2	127,9	1,9	39	122,7	17,9	110	53,8	110,0	Mnat	100
3	453,5	1,6	325	120,9	21,4	23	53,8	108,2	Mnat	99
4	482,1	2,2	28	50,7	32,5	46	39,2	47,7	Bbar	99
5	502,4	1,6	20	120,9	20,5	24	55,8	108,6	Mnat	99
6	550,9	1,5	48	114,9	20,9	23	54,2	112,0	Mnat	99
7	599,9	1,6	48	113,3	23,5	109	57,9	106,2	Mnat	99
8	662,1	1,7	62	107,2	23,9	25	49,0	46,8	Mnat	95
9	725,4	1,7	63	116,6	24,4	106	59,1	106,8	Mnat	100
10	772,9	1,8	47	109,3	17,8	21	48,5	92,6	Mnat	98
11	847,5	1,9	74	113,8	22,4	104	53,8	108,2	Mnat	100
12	905,1	1,5	57	110,2	22,4	24	55,2	97,1	Mnat	98
13	966,8	1,9	61	111,1	15,9	104	47,8	102,0	Mnat	99
14	1025,5	1,8	58	117,6	20,9	108	54,3	108,1	Mnat	100
15	1077,8	1,6	52	116,9	24,5	108	60,9	107,7	Mnat	100
16	1134,6	1,7	56	118,5	22,1	109	58,8	109,4	Mnat	100
17	1181,5	1,7	46	115,6	11,7	16	43,1	22,6	Spec.	0
18	1356,3	1,6	174	120,1	9,0	12	37,5	19,8	Spec.	0
19	1805,9	1,7	449	65,0	22,4	49	39,0	48,9	Spec.	0
20	1859,2	1,7	53	69,2	24,0	47	40,1	48,2	Mnat	78

## 11. CSV documents

**Pro:** When measuring and extracting calls csv files can be generated for use with batIdent and automated identification. These can be opened and displayed in bcAnalyze. A simple graph showing a scatterplot of measurements can be opened as well.

## 12. Preferences

The behavior of bcAnalyze can be adjusted via application wide preferences. These preferences are separated into various topics.

### 12.1. General

The general settings pane holds common settings for bcAnalyze. You can choose a standard channel for display when opening stereo recordings. The delay factor of time-delayed recordings is set here as well. If calls are available for a recording the call list display can be adjusted here as well.

**Pro:** Choose if the first call is automatically selected and zoomed to upon opening a file or if the call list should be displayed automatically if calls are available. If you create a new document you can choose how it should be pre-filled. When exporting measurements to the clipboard a decimal format can be specified.

### 12.2. Display

For the main window of each document the waveform display and the overview sonogram can be adapted.

**Pro:** The pro version allows to make inactive documents semi-transparent and larger fonts to be used for displaying measurements and values.

### 12.3. Call finder

Call finder settings are adjusted here.

### 12.4. Sonogram

Settings for the measurement sonogram (displayed in the extra window) can be adjusted. This includes window size, type and overlap as well as the color scheme.

**Pro:** In addition to the application supplied color schemes you can choose a .tif file with your own color blend. The file must be 1 pixel high and 256 pixels wide.

